Date Received

Date Opened

Kartoffeldextrose-Agar

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8X00131 (1006)

Difco™ Potato Dextrose Agar

Base for Culturing Yeasts and Molds from Food and Dairy Products

Meets USP/EP/JP performance specifications, where applicable.

Directions: Suspend 39 g of the powder in 1 L of purified water. Mix thoroughly. Heat with frequent agitation and boil for 1 minute to completely dissolve the powder. Autoclave at 121°C for 15 minutes. To alter the reaction to pH 3.5, cool the base to 45-50°C and aseptically add an appropriate amount of sterile 10% tartaric acid to each liter of medium. Mix well. Do not reheat the medium. Test samples of the finished product for performance using stable, typical control cultures.

Approximate Formula* Per Liter

Potato Starch (from infusion)** 4.0 g
Dextrose 20.0 g
Agar 15.0 g

*Adjusted and/or supplemented as required to meet performance criteria.

**Approximates 200 g of infusion from potatoes.

For Laboratory Use • Final pH 5.6 ± 0.2

Hygroscopic • Keep container tightly closed.

BD and BD Logo are trademarks of Becton, Dickinson and Company. Difco is a trademark of Difco Laboratories, subsidiary of Becton, Dickinson and Company.

Becton, Dickinson and Company
Sparks, MD 21152 USA • 38800 Le Pont de Claix, France

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Formula approssimativa* per litro

Amido di patata (infuso)** 4,0 g
Destrosio 20,0 g
Agar 15,0 g

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