# Extendable Pole

English

## 3M™ Extendable Pole

**REF** XSL008

**Description:** Extendable aluminum pole with Sponge-Stick holder, 2.56 m (8 ft 5 in)

**Applications:** For Laboratory Use Only.

### Instructions for Use

**(Additional requirement: REF SSL100, SSL10BPW, SSL10LETG, SSL10NB or SSL10NBG)**

- 1) Loosen the screw on the stick holder section of the pole.
- 2) Shake stick to end of bag.
- 3) Tear bag open.
- 4) Aseptically, push stick out of bag until thumb-stop line is exposed. Remove Sponge-Stick from bag.
- 5) Insert stick into holder. Secure stick with screw.
- 6) Turn each pole section in opposite direction to release lock. Pull out to preferred length. Turn pole section in opposite direction to lock.
- 7) Aseptically swab across the entire sampling surface. (See Caution below.)
- 8) Turn sponge over. Change direction 90°. Aseptically swab the same sampling surface.
- 9) Turn each pole section in opposite direction to release lock. Push in to collapse. Turn pole section in opposite direction to lock.
- 10) Loosen screw to remove Sponge-Stick from holder.
- 11) Aseptically insert sponge head into bag.
- 12) Hold sponge in place inside bag. Bend stick to break. Allow sponge to drop in bag. Discard stick.
- 13) Fold bag to close.
- 14) Fold ends of blue wire inward.
- 15) Retain pole for future use.

**CAUTION:** DO NOT USE NEAR ELECTRICAL WIRES.

**Disposal:** Follow established laboratory practices.


### 3M

St. Paul, MN, 55144-1000 USA

[www.3M.com/microbiology](http://www.3M.com/microbiology)

Made in U.S.A.

3M is a trademark of 3M and used under license in Canada. ©3M 2008. All rights reserved.  
34-8700-5597-6

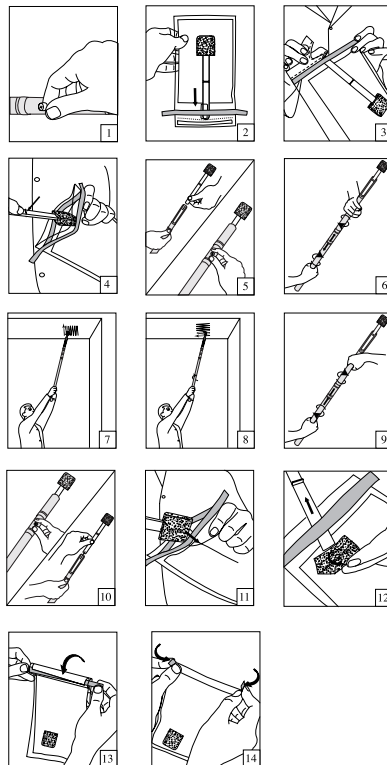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Trademark 3M es una marca registrada de 3M y se usa bajo licencia en Canadá. ©3M 2008.