

3M Food Safety Environmental Listeria Monitoring Solution



Zero In On *Listeria*



3M™ Petrifilm™
Environmental Listeria Plate

3M™ Mini Flip-Top Vial w/10mL
Buffered Peptone Water Broth

3M™ Quick Swab


Monitor environmental *Listeria* levels with consistent, frequent testing to identify hot spots in your plant and track your progress. Cost effectively test more often in more locations. Zero in on the source of *Listeria* with a complete solution from 3M Food Safety.

- Quantitative results allow you to identify hot spots in the plant
- Rapid results confirmed in 27–31 hours from sample collection
- Safer method—no enrichment, no culture transfers—less risk of cross-contamination
- No specialized equipment needed
- Easy-to-use solution to make faster decisions, track progress against *Listeria* over time, and react to problem areas







Fast, Easy and More Informative

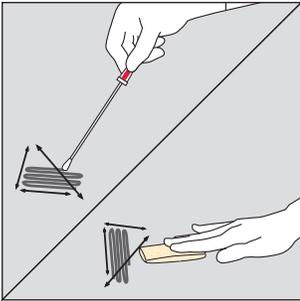


1. Sample using swab or sponge.

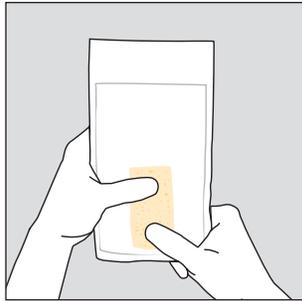

2. Add buffered peptone water. Wait one hour.

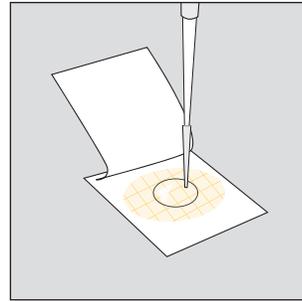

3. Inoculate and incubate for 26–30 hours.

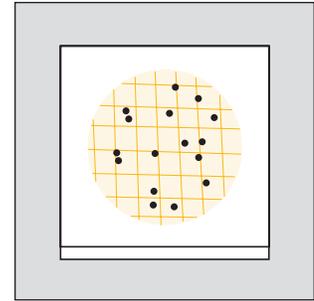

4. Read and interpret.

End-to-end solution that works together for consistent, reliable results

Product Solutions

Microbial Testing—For detection of *Listeria* in environmental samples

Product	Catalog Number	Quantity
3M™ Petrifilm™ Environmental Listeria Plate	6447	50 plates
	6448	200 plates

Sample Collection

Product	Catalog Number	Quantity
3M™ Quick Swab	6432	50 swabs
	6433	250 swabs
3M™ Swab-Sampler w/1mL Letheen Broth	RS9601LET	100 swabs
3M™ Swab-Sampler w/4mL Letheen Broth	RS9604LET	100 swabs
3M™ Swab-Sampler w/5mL Letheen Broth	RS9605LET	100 swabs
3M™ Swab-Sampler w/10mL Letheen Broth	RS96010LET	100 swabs
3M™ Enviro Swab	ENWSWB25	25 swabs
3M™ Sponge-Stick, 1.5" x 3" Sponge with 10mL Letheen Broth	SSL10LET	100 sponges (5 bags of 20)
3M™ Hydrated-Sponge, 1.5" x 3" Sponge with 10mL Letheen Broth	HS10LET	100 sponges (5 bags of 20)

Repair and Recovery—For repair of *Listeria* in environmental samples

Product	Catalog Number	Quantity
3M™ Mini Flip-Top Vial w/10mL Buffered Peptone Water Broth	BPPFV10BPW	100 vials
Buffered Peptone Water for repair of <i>Listeria</i> in environmental samples, 200mL	6449	20 vials



3M Food Safety

3M Center
Building 275-5W-05
St. Paul, MN 55144-1000
USA
1-800-328-6553
www.3M.com/foodsafety

Please recycle. Printed in U.S.A.
© 3M 2011. All rights reserved.
70-2009-9498-9

3M and Petrifilm are trademarks of 3M.

For more information visit us at
www.3M.com/foodsafety or call 1-800-328-6553